# Product Description

# **SALSA® Binning DNA SD078-S01**

#### **Version S01**

#### Catalogue number

• SD078: SALSA Binning DNA, 6 reactions

#### **Precautions and warnings**

For professional use only. Always consult the most recent product description AND the corresponding probemix product description AND the MLPA General Protocol before use: www.mrcholland.com. Binning DNA is not known to contain any harmful agents.

#### Safety data sheet

Based on the concentrations present, none of the ingredients are hazardous as defined by the Hazard Communication Standard. A Safety Data Sheet (SDS) is not required for these products: none of the preparations contain dangerous substances (as per Regulation (EC) No 1272/2008 [EU-GHS/CLP] and amendments) at concentrations requiring distribution of an SDS (as per Regulation (EC) No 1272/2008 [EU-GHS/CLP] and 1907/2006 [REACH] and amendments). If spills occur, clean with water and follow appropriate site procedures.

#### **General information**

The SALSA Binning DNA SD078 is a research use only (RUO) reagent to be used in combination with SALSA MLPA Probemix P190-D1 CHEK2, a SALSA MLPA Reagent Kit and Coffalyser.Net™ analysis software for the processes of linking all probe signals to their identity by use of the probe lengths. SD078 contains the targets of all probes included in the above-listed probemix, including the mutation-specific probe target CHEK2 c.1100delC.

Binning DNA should never be used as a reference sample in the MLPA data analysis. Neither should it be used in quantification of mutation signals.

#### **Experimental set up**

MLPA reactions for binning purposes should be performed with  $5\,\mu$ l of Binning DNA. Inclusion of one reaction with SALSA Binning DNA SD078 in the initial MLPA experiment is essential as it can aid in data binning of the peak pattern when using Coffalyser.Net software. Furthermore, Binning DNA should be included in the experiment whenever changes have been applied to the set-up of the capillary electrophoresis device (e.g. when a different polymer type is used).

#### **Data analysis**

Coffalyser.Net software should be used for analysis of MLPA experiments. When performing the fragment analysis step in Coffalyser.Net, select SD078 in the *bin smpl* –column. By selecting the SD078 sample as your binning sample, probes will be correctly identified in the peak pattern across all samples. Coffalyser.Net software is freely downloadable at <a href="https://www.mrcholland.com">www.mrcholland.com</a>.

### **Binning DNA content**

SD078 consists of a mixture of female genomic DNA from healthy individuals and a titrated amount of synthetic DNA that contains a partial sequence of the *CHEK2* gene. This partial sequence includes one mutation that will be detected by the mutation-specific probe present in the above-listed probemix. See Table 1 and the corresponding probemix product description for more details on the mutation-specific probe target present. The indicated mutation-specific probe will generate a signal on SD078.

Please note that the synthetic DNA also contains the target sequence of the 105 nt chromosome Y specific control fragment. As a result, the 100 and 105 nt control fragments indicate the presence of two copies chromosome X and one copy chromosome Y.

SALSA Binning DNA SD078



## Table 1. Mutation-specific probe target in Binning DNA SD078-S01

Probemix	Gene/Exon	Probe length (nt)	Probe ID	Probemix version	Details
P190	CHEK2 exon 11	313	22034-SP0468-L31261	D1	c.1100delC

**Note:** Please consult the corresponding probemix product description for more information about exon numbering, mutation nomenclature and gene transcripts used.

More information: www.mrcholland.com; www.mrcholland.eu				
***	MRC Holland bv; Willem Schoutenstraat 1			
	1057 DL, Amsterdam, The Netherlands			
E-mail	info@mrcholland.com (information & technical questions)			
	order@mrcholland.com (orders)			
Phone	+31 888 657 200			

IVD	EUROPE* (until lot 1021)	CE
RUO	EUROPE* (as of lot 0325) ALL OTHER COUNTRIES	

<sup>\*</sup>comprising EU (candidate) member states and members of the European Free Trade Association (EFTA), and the UK. The product is for RUO in all other European countries.

#### Implemented changes in the product description

Version S01-05 - 01 May 2025 (03)

- Product description updated as SD078 is for research use only from lot 0325 onwards.
- CHEK2 mutation nomenclature modified from "1100delC" to "c.1100delC".

Version S01-04 - 22 October 2021 (03)

- Product description adapted to a new template.
- Intended purpose updated.

Version S01-03 - 30 October 2020 (02)

- Product description adapted to a new template.
- Intended purpose updated.

SALSA Binning DNA SD078 Page 2 of 2